# lechnotes

Highlighting technical features for **Thermo Scientific Vacutherm Vacuum Ovens** 

#### **Key Features**

- Three chamber sizes available: 25L, 53L, 128L
- Outstanding capabilities for fast and gentle drying
- Customizable options available to fit your application needs

## Why choose a Vacutherm vacuum oven for your lab?

Dry up to six times faster than conventional units with a Thermo Scientific<sup>™</sup> Vacutherm<sup>™</sup> vacuum heating and drying oven.\* Minimize operation time with outstanding heat transfer and fast heat up.\*\*



#### MODELS

- Jacket-heated models, available with Kelvitron<sup>™</sup> microprocessor control, operate up to 200°C
- Shelf heated models operate up to 300°C (400°C option offered), available with Digicon multi-channel controls which allow you to manage each shelf temperature separately

#### **TYPICAL APPLICATIONS**

#### Food industry

- Determination of the solid content of foods. Samples are typically dried under vacuum at around 60°C
- ~ Our Solution: Vacutherm VT6060 M or VT6130 M oven with standard pump and trap Pharmaceutical industry
- Drying additives for pharmaceuticals. Samples are dried with pressure at 10-1mbar or forming gas at 60-80°C.
- ~ Our Solution: Vacutherm VT6060 P or VT6130 P oven with an oil-free pump
- Gentle drying of raw materials and determination of solid content. Samples are dried under vacuum at 60-80°C.
- ~ Our Solution: Vacutherm VT6060 M or VT6130 M oven with standard pump and trap
- \* Based on internal testing
- \*\* Based on previous models

| Specifications and Ordering Information |           |  |                       |   |   |                                |  |                                |                        |                    |
|---|-----------|--|-----------------------|---|---|--------------------------------|--|--------------------------------|------------------------|--------------------|
| Cat.No.                                 | Model     | Туре   | Capacity              | Interior Dimensions<br>W x H x D              | Exterior Dimensions <sup>1</sup><br>W x H x D | Rated<br>Temp. <sup>2, 3</sup> | Spatial temp.<br>deviation <sup>2, 4</sup> | Heat Output<br>(at max. temp.) | Electrial <sup>5</sup> | Weight             |
| 51014550                                | VT 6025   | Jacket-heated<br>no inert gas<br>connection  | 25L<br>(0.88 cu.ft.)  | 300 x 275 x 307mm<br>(11.8 x 10.8 x 12.1in.)  | 480 x 600 x 450 mm<br>(18.9 x 23.6 x 17.7in.) | 200°C                          | ±4 (@200°C)                                | 340 Wh/h                       | 230V 50/60Hz,<br>1.3kW | 58kg<br>(128 lb.)  |
| 51014552                                | VT 6025   | Jacket-heated                                | 25L<br>(0.88 cu.ft.)  | 300 x 275 x 307mm<br>(11.8 x 10.8 x 12.1in.)  | 480 x 600 x 450 mm<br>(18.9 x 23.6 x 17.7in.) | 200°C                          | ±4 (@200°C)                                | 340 Wh/h                       | 230V 50/60Hz,<br>1.3kW | 58kg<br>(128 lb.)  |
| 51014553                                | VT 6025   | Jacket-heated6                               | 25L<br>(0.88 cu.ft.)  | 300 x 275 x 307mm<br>(11.8 x 10.8 x 12.1in.)  | 480 x 600 x 450 mm<br>(18.9 x 23.6 x 17.7in.) | 200°C                          | ±4 (@200°C)                                | 340 Wh/h                       | 230V 50/60Hz,<br>1.3kW | 58kg<br>(128 lb.)  |
| 51014554                                | VT 6025   | Jacket-heated<br>with Digicon S <sup>7</sup> | 25L<br>(0.88 cu.ft.)  | 300 x 275 x 307mm<br>(11.8 x 10.8 x 12.1in.)  | 480 x 600 x 450 mm<br>(18.9 x 23.6 x 17.7in.) | 200°C                          | ±4 (@200°C)                                | 340 Wh/h                       | 230V 50/60Hz,<br>1.3kW | 58kg<br>(128 lb.)  |
| 51014539                                | VT 6060 M | Jacket-heated                                | 53L<br>(1.87 cu.ft.)  | 415 x 345 x 371 mm<br>(16.3 x 13.6 x 14.6in.) | 744 x 576 x 570 mm<br>(29.3 x 22.7 x 22.4in.) | 200°C                          | ±4 (@200°C)                                | 540Wh/h                        | 230V 50/60Hz,<br>1.7kW | 82kg<br>(181 lb.)  |
| 51014542                                | VT 6060 P | Shelf-heated <sup>8</sup>                    | 53L<br>(1.87 cu.ft.)  | 415 x 345 x 371 mm<br>(16.3 x 13.6 x 14.6in.) | 744 x 576 x 570 mm<br>(29.3 x 22.7 x 22.4in.) | 300/400°C                      | ±3 (@200°C)<br>±7 (@300°C)<br>±9 (@400°C)  | 450 Wh/h                       | 230V 50/60Hz,<br>1.6kW | 90kg<br>(198 lb.)  |
| 51014541                                | VT 6130 M | Jacket-heated                                | 128L<br>(4.52 cu.ft.) | 495 x 489 x 529<br>(19.5 x 19.3 x 20.8in.)    | 895 x 720 x 750 mm<br>(35.2 x 28.3 x 29.5in.) | 200°C                          | ±7 (@200°C)                                | 870 Wh/h                       | 230V 50/60Hz,<br>2.2kW | 153kg<br>(337 lb.) |
| 51014543                                | VT 6130 P | Shelf-heated <sup>8</sup>                    | 128L<br>(4.52 cu.ft.) | 495 x 489 x 529 mm<br>(19.5 x 19.3 x 20.8in.) | 895 x 720 x 750 mm<br>(35.2 x 28.3 x 29.5in.) | 300/400°C                      | ±4 (@200°C)<br>±7 (@300°C)<br>±9 (@400°C)  | 880 Wh/h                       | 230V 50/60Hz,<br>3.0kW | 164kg<br>(362 lb.) |

1 Depth without handle and wall distance. Dimensions: handle 70 mm, wall distance 80 mm 2 The values stated apply to the empty oven and vacuum operation in acc. with DIN 12880, part 2

3 Control range M: electronic controller from TA + 15°C; control range P: electronic controller from TA +10°C; TA = Ambient temperature

4 Measured on the shelves

5 Other voltages on request 6 With digital pressure display and recorder output

7 Digicon S: analog temperature controller

8 Comes standard with Digicon multi-channel controller and rated temperature of 300°C; for 400°C: order option in addition, see next page



### Find the right Vacutherm vacuum oven for your application needs



- Gentle drying of heat sensitive materials
- Significantly reduced drying times when compared to convectional oven without vacuum
- Drying of intricately designed parts
- Elimination of oxidation associated with heat treatment
- Targeted recovery of degassed products

|  | ROTARY<br>Vane Pumps |                 |                 |                 |                 |  |
|--|----------------------|-----------------|-----------------|-----------------|-----------------|--|
| Vacuum<br>Pumps  |                      | MD 12H          | HMD 4C          | HMZ 2C          | RE 9H           |  |
| Cat. No.   |                      | 50040078        | 50028362        | 50028364        | 50028361        |  |
| Rated pumping<br>capacity m <sub>3</sub> /h<br>(Pneurop) |                      | 9.6 3.0         |                 | 1.7             | 8.6             |  |
| Total final<br>pressure without<br>gas ballast           | mbar                 | 2               | 2               | 10              | 0.1             |  |
| Total final<br>pressure with<br>gas ballast              | mbar                 | <10             | <10             | <20             | 6 x 10-1        |  |
| Water vapor<br>compatibility                             | mbar                 | _               | _               | _               | 40              |  |
| Capacity of trap   | ml                   | 725             | 500             | 500             | 500             |  |
| Weight   | kg                   | 24              | 18              | 14              | 21.6            |  |
| Dimensions with trap (w/h/d)                             | mm                   | 240 x 300 x 610 | 241 x 500 x 338 | 241 x 326 x 336 | 260 x 232 x 460 |  |
| Vacucenter<br>with pump                                  |                      | MD 12H          | HMD 4C          |                 | RE 9H           |  |
| Cat. No.*  |                      | 50044332        | 50044333        | -               | 50044334        |  |
| Voltage  | V3N~,<br>50/60 Hz    | 400             | 400             | _               | 400             |  |
| Max. rated<br>power                                      | kW                   | 3.37            | 3.2             | -               | 3.37            |  |
| External<br>dimensions<br>(W x H x D)                    | mm                   | 895 x 650 x 900 | 895 x 650 x 900 | -               | 895 x 650 x 900 |  |
| Total weight   | kg                   | 88              | 80              | -               | 86              |  |
| Protection class   |                      | 1               | 1               | -               | 1               |  |
| Protection type  |                      | IP 20           | IP 20           | -               | IP 20           |  |

\*In combination with VT 6060 or VT 6130

| ACCESSORI  | ES   |  |  |  |  |
|------------|--|--|--|--|--|
| Cat.No.    | Description  |  |  |  |  |
| 50028403   | Additional shelf for VT6025 (incl. shelf supports)   |  |  |  |  |
| 50043975   | Additional shelf for VT6060 M (incl. shelf supports)   |  |  |  |  |
| 50043976   | Additional shelf for VT6130 M (incl. shelf supports)   |  |  |  |  |
| 50048621   | Stainless steel tray for VT6060  |  |  |  |  |
| 50048619   | Stainless steel tray for VT6130  |  |  |  |  |
| 50048620   | Aluminum tray for VT6060 M/P   |  |  |  |  |
| 50048618   | Aluminum tray for VT6130 M/P   |  |  |  |  |
| 50046860   | Stainless steel vacuum connection kit for tubing, 10mm diameter (all models)                                     |  |  |  |  |
| 50029890   | Support frame, 780 mm height, for VT6060 M/P   |  |  |  |  |
| 50029597   | Support frame, 780 mm height, for VT6130 M/P   |  |  |  |  |
| FACTORY IN | ISTALLED OPTIONS   |  |  |  |  |
| Cat.No.    | Options for Vacutherm 6060 and 6130 M/P  |  |  |  |  |
| 51900284   | RS 232 computer interface for Kelvitron controller   |  |  |  |  |
| 51900069   | Digital pressure display   |  |  |  |  |
| 51900193   | Pressure controller with solenoid valve  |  |  |  |  |
| 51900071   | VITON door gasket for VT 6060 M  |  |  |  |  |
| 51900072   | VITON door gasket for VT 6130 M  |  |  |  |  |
| 51900205   | Daily program timer, mechanical  |  |  |  |  |
| 51900073   | 24 hour timer (20 h at 60 Hz)  |  |  |  |  |
| 51900008   | Weekly program timer, mechanical   |  |  |  |  |
| 51900161   | Weekly program timer, digital  |  |  |  |  |
| 51900074   | Digital sample temperature display with flexible pt 100 temperature sensor and socket for external data recorder |  |  |  |  |
| 51900209   | Thermicon P temperature program controller (M-models only)   |  |  |  |  |
| 51900297   | Eurotherm 2404/P4 temperature controller with RS232 interface (M-models only)                                    |  |  |  |  |
| 51900298   | Eurotherm 2404/P4 temperature controller with RS 422/485 interface (M-models only)                               |  |  |  |  |
| 51900081   | Central monitoring connection  |  |  |  |  |
| Cat.No.    | Digicon <sup>®</sup> S temperature controller with analog interface 0–20 mA/0–10 V                               |  |  |  |  |
| 51900194   | (200 °C): VT 6060 M and VT 6130 M  |  |  |  |  |
| 51900195   | (300 °C): VT 6060 P and VT 6130 P  |  |  |  |  |
| 51900196   | (400 °C): VT 6060 P  |  |  |  |  |
| 51900197   | (400 °C): VT 6130 P  |  |  |  |  |
| Cat.No.    | Rated temperature of 400 °C, (door without inspection window)  |  |  |  |  |
| 51900079   | for VT 6060 P  |  |  |  |  |
| 51900080   | for VT 6130 P  |  |  |  |  |
| Cat. No.   | Vacuum chamber with rounded corners at the rear  |  |  |  |  |
| 51900075   | for VT 6060 M  |  |  |  |  |
| 51900077   | for VT 6060 P  |  |  |  |  |
| 51900076   | for VT 6130 M  |  |  |  |  |
| 51900078   | for VT 6130 P  |  |  |  |  |
| Cat. No.   | Options for Vacutherm 6025   |  |  |  |  |
| 51900326   | Precision control valve for standard model (order nos. 51014550 and 51014551 only)                               |  |  |  |  |
| 51900284   | RS232 computer interface   |  |  |  |  |

#### thermoscientific.com/vacuumovens

© 2014 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Australia +61 39757 4300 Austria +43 1 801 40 0 Belgium +32 53 73 42 41 China +800 810 5118 or +400 650 5118 France +33 2 2803 2180 Germany national toll free 0800 1 536 376 Germany international +49 6184 90 6000 India toll free 1800 22 8374 India +91 22 6716 2200 Italy +39 02 95059 552 Japan +81 3 5826 1616 Netherlands +31 76 579 55 55 New Zealand +64 9 980 6700 Nordic/Baltic/CIS countries +358 9 3291 0200 Russia +7 812 703 42 15 Spain/Portugal +34 93 223 09 18 Switzerland +41 44 454 12 22 UK/Ireland +44 870 609 9203 USA/Canada +1 866 984 3766

**Other Asian countries** +852 2885 4613 **Countries not listed** +49 6184 90 6000

